

Mouse Anti-Type II Collagen IgG Antibody Subtype ELISA Kits with TMB

Catalog # 20311T-20315T, 20321T-20325T, 20331T-20335T, 20351T-20355T, 20361T-20365T

For Research Use Only - Not Human or Therapeutic Use

PRODUCT SPECIFICATIONS

DESCRIPTION:	Assay kit to quantify mouse anti-collagen antibodies
FORMAT:	Pre-coated 96-well ELISA Plate with removeable strips
ASSAY TYPE:	Indirect ELISA
ASSAY TIME:	4.25 hours
STANDARD RANGE:	IgG1, IgG2a, IgG2b, and IgG2c : 10-0.16 ng/ml IgG3 : 100-1.6 ng/ml
NUMBER OF SAMPLES:	Up to 39 (duplicate) samples/standard plate (will vary for custom kits)
SAMPLE TYPES:	Serum and Plasma
RECOMMENDED SAMPLE DILUTIONS:	1:1000 (at least)
CHROMOGEN:	TMB (read at 450 nm)
STORAGE:	-20°C
VALIDATION DATA:	N/A
NOTES:	

Mouse Anti-Type II Collagen IgG Antibody Subtype ELISA Kits with TMB

Catalog # 20311T-20315T, 20321T-20325T, 20331T-20335T, 20351T-20355T, 20361T-20365T

For Research Use Only - Not Human or Therapeutic Use

INTRODUCTION

Autoantibodies to type II collagen play a primary role in the collagen induced arthritis (CIA) model. However, autoantibodies are not always capable of inducing arthritis due to the inability to activate a complement, the first critical step in the activation of inflammatory cascades. In CIA susceptible DBA mice, complement fixable IgG2a and IgG2b subtype autoantibody levels, as well as their epitope specificity, are pertinent in the induction of arthritis. In order to adequately study the pathological roles of antibodies to type II collagen, it is important to determine individual IgG subtype autoantibody levels to mouse type II collagen as well as the heterologous type II collagen used for immunization. On the other hand, C57BL/6 mice, a popular strain for developing transgenic mice, produce IgG2c antibodies instead of IgG2a antibodies due to its MHC class I haplotype (H-2b). Moreover, these mice can only develop CIA when immunized with chick type II collagen. Therefore, serum IgG2c autoantibody levels against mouse type II collagen and chick type II collagen should be evaluated when using C57BL/6 mice for the CIA model.

Chondrex, Inc. provides Mouse Anti-Collagen IgG Antibody Subtype (IgG1, IgG2a, IgG2b, IgG2c, and IgG3) ELISA Kits for further analysis of antibodies in mouse serum from the mouse CIA model. Depending on the IgG subtypes and epitopes, the antibody-antigen affinity varies significantly among individual antibodies and serum samples from individual animals. Therefore, the total IgG concentration calculated as the sum of individual IgG subtypes might not perfectly match the total IgG levels determined by Chondrex, Inc's Mouse Anti-Collagen IgG ELISA Kit. These kits can also be used to determine antibody levels to type I collagen (chick, bovine, porcine, human, and mouse) upon request. Please contact Chondrex, Inc. at support@chondrex.com for more information on custom type I collagen-coated plates.

KIT COMPONENTS

	Item	Quantity	Amount	Storage
Standard Antibody	IgG1 (20306) -10 ng	1 vial	Lyophilized	-20°C
	IgG2a (20307) - 10 ng			
	IgG2b (20308) -10 ng			
	IgG2c (20309) -10 ng			
	IgG3 (20305) - 100 ng			
Peroxidase-Conjugated Anti-Mouse Secondary Antibody	IgG1 (203011)	2 vials	50 µl	-20°C
	IgG2a (203012)			
	IgG2b (203013)			
	IgG2c (203010)			
	IgG3 (203017)			
	Solution A - Blocking Buffer (2071)	1 bottle	10 ml	-20°C
	Solution B - Sample/Standard Dilution Buffer (2072)	1 bottle	50 ml	-20°C
	Solution C - Secondary Antibody Dilution Buffer (2073)	1 bottle	20 ml	-20°C
	TMB (90023)	2 vials	0.2 ml	-20°C
	Chromogen Dilution Buffer (90022)	1 bottle	20 ml	-20°C
	Stop Solution - 2N Sulfuric Acid (9016)	1 bottle	10 ml	-20°C
	Wash Buffer, 20X (9005)	1 bottle	50 ml	-20°C
	Type I or Type II Collagen Coated 8-Well Strips	10 each	8-well strips	-20°C
	Reference Standard Strips (two strips per run)	4 each	8-well strips	-20°C


PLATE COATING AND SETUP

Species (Type II Collagen Color Coding)	IgG1 Catalog #	IgG2a Catalog #	IgG2b Catalog #	IgG2c Catalog #	IgG3 Catalog #
Chick (Yellow)	20311T	20312T	20313T	20315T	20314T
Bovine (Green)	20321T	20322T	20323T	20325T	20324T
Porcine (Pink)	20331T	20332T	20333T	20335T	20334T
Human (Blue)	20351T	20352T	20353T	20355T	20354T
Mouse (Orange)	20361T	20362T	20363T	20365T	20364T
Standard (Red)	---	---	---	---	---


Standard ELISA Kit with One Species of Type II Collagen

A standard ELISA kit consisting of ten 8-well strips coated with one species of type II collagen and two 8-well strips for reference standards. "B" represents blank wells to determine background values caused by the secondary antibody. Standards and samples (1-39) are run in duplicate.

Note: custom coating of plates is available upon request.

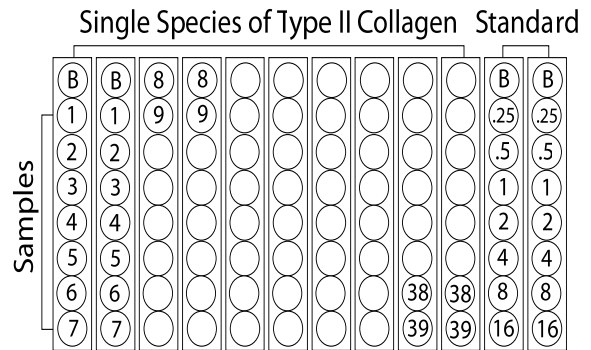

ASSAY OUTLINE

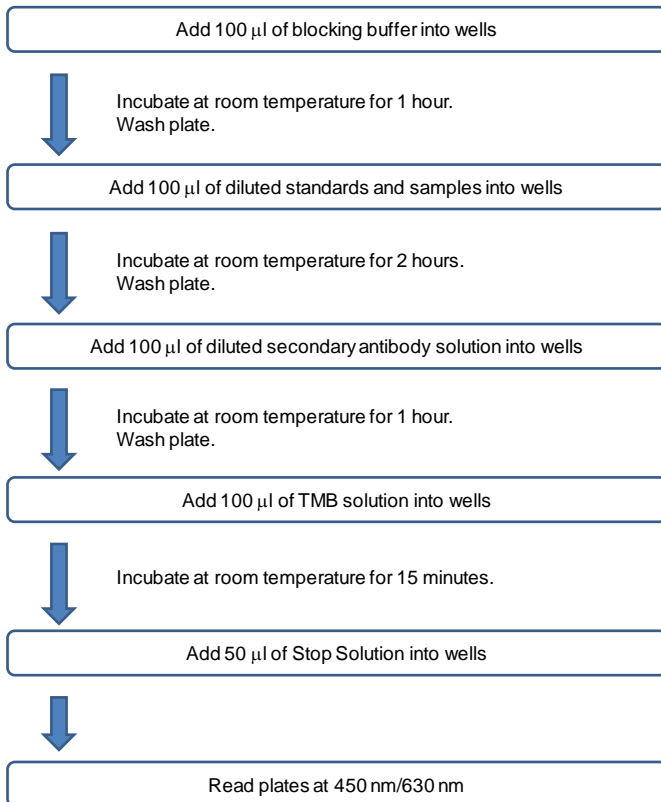

NOTES BEFORE USING ASSAY

NOTE 1: It is recommended that the standard and samples be run in duplicate.

NOTE 2: Warm up all buffers to room temperature before use.

NOTE 3: Crystals may form in Wash Buffer, 20X when stored at cold temperatures. If crystals have formed, warm the wash buffer by placing the bottle in warm water until crystals are completely dissolved.

NOTE 4: Measure exact volume of buffers using a serological pipet, as extra buffer is provided.


NOTE 5: Cover the plate with plastic wrap or a plate sealer after each step to prevent evaporation from the outside wells of the plate.

NOTE 6: For partial reagent use, please see the assay protocol's corresponding step for the appropriate dilution ratio. For example, if the protocol dilutes 50 µl of a stock solution in 10 ml of buffer for 12 strips, then for 6 strips, dilute 25 µl of the stock solution in 5 ml of buffer. Partially used stock reagents may be kept in their original vials and stored at -20°C for use in a future assay.

NOTE 7: This kit contains animal components from non-infectious animals and should be treated as potential biohazards in use and for disposal.

ASSAY PROCEDURE

- Add Blocking Buffer:** Add 100 µl of Blocking Buffer (Solution A) to all wells. Incubate for 1 hour at room temperature.
- Prepare Standard Dilutions:** Dissolve one vial of standard with 1 ml of Sample/Standard Dilution Buffer (Solution B) to make the highest standard concentration, labeled "1" in the standard dilution diagram below. Prepare serial dilutions of the 1X standard by mixing 250 µl of the 1X standard with 250 µl of Solution B, labeled "2" in the standard dilution diagram below. Then repeat this procedure five more times to make a total of seven standard stock solutions. The 1X standard may be stored at -20°C for use in a second assay. Chondrex, Inc. recommends making fresh serial dilutions for each assay.

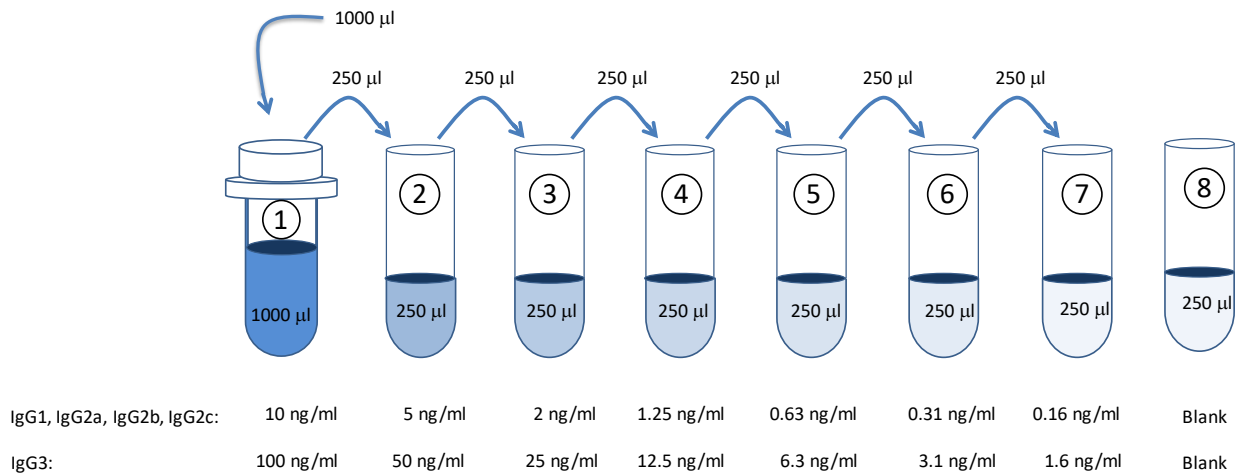

3. **Prepare Sample Dilutions:** Centrifuge serum samples at 10,000 rpm at room temperature for 3 minutes to remove insoluble materials and lipids. Dilute samples 1:1000 or more with Solution B. For example, dilute 10 μ l of sample with 0.99 ml of Solution B (1:100). Keep this as a stock solution for future assays. If it is necessary to assay antibodies at a low dilution (less than 1:200) due to low antibody levels, please contact Chondrex, Inc. customer service at support@chondrex.com. In the CIA model, anti-collagen antibody levels may vary from zero (naïve rodents) to 1,000,000 units/ml or higher (arthritic rodents) depending on the immune response, arthritis severity, and time of sample collection. With unknown anti-collagen antibody levels, Chondrex, Inc. recommends testing several dilutions, such as 1:1000, 1:10,000, and 1:100,000, in order to find the appropriate dilution for the samples.
4. **Dilute Wash Buffer:** Dilute 50 ml of 20X wash buffer in 950 ml of distilled water (1X wash buffer). Wash the plate with 1X wash buffer at least 3 times using a wash bottle with manifold or an automated plate washer. Empty the plate by inverting it and blotting on a paper towel to remove excess liquid. *Do not allow the plate to dry out.*
5. **Add Standards and Samples:** Add 100 μ l of standards, Solution B (blank) and samples to wells in duplicate. Incubate at room temperature for 2 hours.
6. **Wash:** Wash the plate with 1X wash buffer at least 3 times using a wash bottle with manifold or an automated plate washer. Empty the plate by inverting it and blotting on a paper towel to remove excess liquid. *Do not allow the plate to dry out.*
7. **Add Secondary Antibody:** Dilute one vial of secondary antibody in 10 ml Secondary Antibody Dilution Buffer (Solution C). Add 100 μ l of secondary antibody solution to each well and incubate at room temperature for 1 hour.

Strip #	Secondary Antibody (μ l)	Solution C (ml)
2	8	1.7
4	17	3.3
6	25	5.0
8	33	6.6
10	42	8.2
12	50	10.0

8. **Wash:** Wash the plate with 1X wash buffer at least 3 times using a wash bottle with manifold or an automated plate washer. Empty the plate by inverting it and blotting on a paper towel to remove excess liquid. *Do not allow the plate to dry out.*
9. **TMB:** Use new tubes when preparing TMB. Dilute one vial of TMB with 10 ml of Chromogen Dilution Buffer just prior to use. Add 100 μ l of TMB solution to each well immediately after washing the plate. Incubate for 15 minutes at room temperature.

Strip #	TMB (μ l)	Chromogen Dilution Buffer (ml)
2	34	1.7
4	66	3.3
6	100	5.0
8	132	6.6
10	164	8.2
12	200	10.0

10. **Stop:** Add 50 μ l of 2N sulfuric acid (Stop Solution) to each well.
11. **Read Plate:** Read the OD values at 450 nm. If the OD values of samples are greater than the OD values of the highest standard, re-assay the samples at a higher dilution. A 630 nm filter can be used as a reference.


CALCULATION OF ANTIBODY TITERS

1. Average the duplicate OD values for the standards, blanks (B) and test samples.
2. Subtract the blank (B) values from the averaged OD values of the standards and test samples.

NOTE: Individual antigens have unique background values. Therefore, blank wells should be used for each different antigen.

3. Plot the OD values of standards against the ng/ml of antibody standard. Using a log/log plot will linearize the data. Figure 1 shows examples of standard curves for IgG1, IgG2a, IgG2b, IgG2c, and IgG3 subtype antibodies.
4. The ng/ml of antibody in test samples can be calculated using regression analysis.

Figure 1 - Typical Standard Curves for the Mouse Anti-Type II Collagen IgG1, IgG2a, IgG2b, IgG2c, and IgG3 Antibody Subtype ELISA Kits.

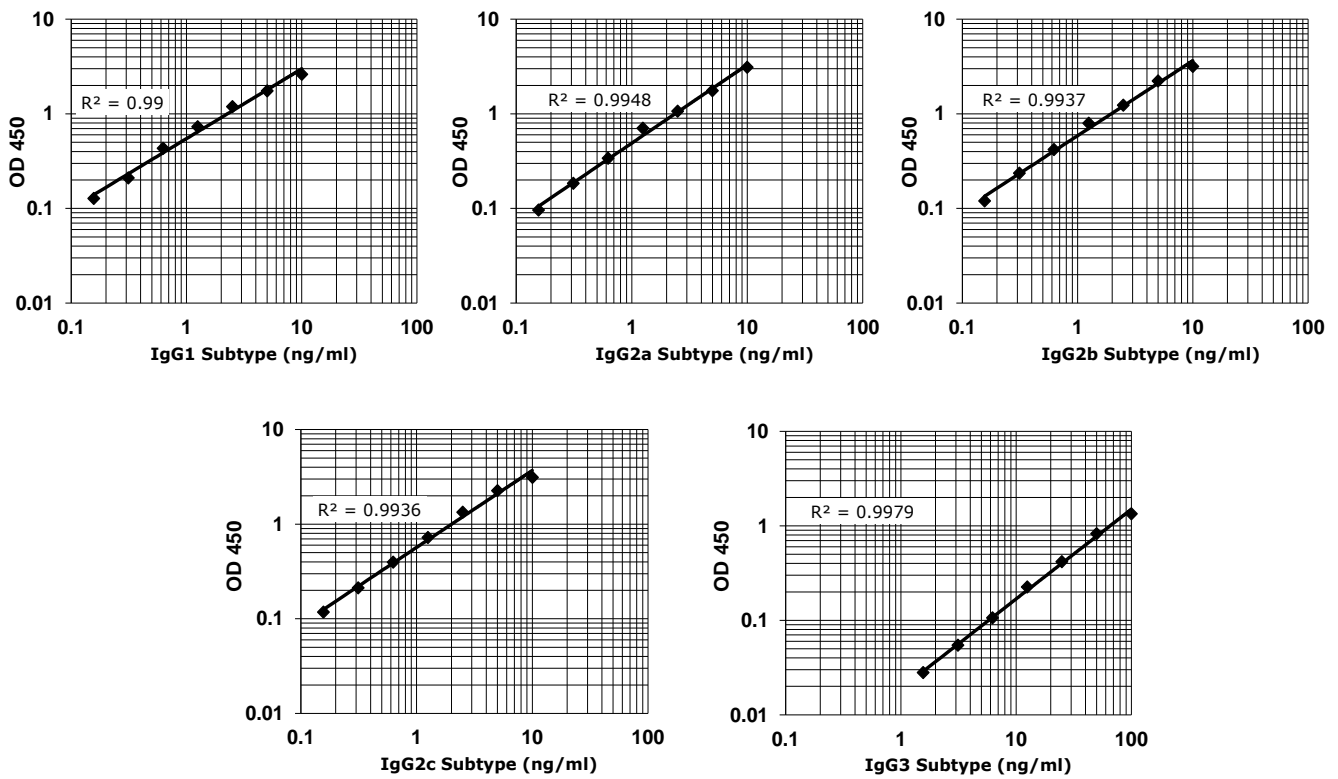

TROUBLESHOOTING

For frequently asked questions about assays and ELISAs, please see Chondrex, Inc.'s [ELISA FAQ](#) for more information.