

Mouse Anti-Glutenin Antibody Subtype/Subclass ELISA Kits

Catalog # 3092, 3093, 3094, 3095, and 3096

For Research Use Only - Not Human or Therapeutic Use

PRODUCT SPECIFICATIONS

DESCRIPTION:	ELISA kit to quantify mouse anti-glutenin antibodies
FORMAT:	96-well ELISA Plate with removeable strips
ASSAY TYPE:	Indirect ELISA
ASSAY TIME:	3.5 hours
STANDARD RANGE:	3092 (IgG) : 100 – 1.6 ng/ml 3093 (IgG1) : 100 – 1.6 ng/ml 3094 (IgG2a) : 100 – 1.6 ng/ml 3095 (IgG2b) : 200 – 3.2 ng/ml 3096 (IgM) : 100 – 1.6 ng/ml
NUMBER OF SAMPLES:	Up to 40 (duplicate) samples/plate
SAMPLE TYPES:	Serum & Plasma
RECOMMENDED SAMPLE DILUTIONS:	1:100 (at least): Depending on immunization protocols
CHROMOGEN:	TMB (read at 450 nm)
STORAGE:	-20°C for 12 months
VALIDATION DATA:	3092: Intra-Assay (2.6-3.1%)/Inter-Assay (3.4-5.3%)/Spiking Test (97-108%) 3093: Intra-Assay (0.9-3.5%)/Inter-Assay (3.7-10%)/Spiking Test (102-108%) 3094: Intra-Assay (6.5-9.1%)/Inter-Assay (1.9-5.9%)/Spiking Test (98-103%) 3095: Intra-Assay (4.8-9.4%)/Inter-Assay (4.5-8.4%)/Spiking Test (97-105%) 3096: Intra-Assay (2.3-7.6%)/Inter-Assay (0.5-5.6%)/Spiking Test (98-109%)
NOTES:	N/A

Mouse Anti-Glutenin Antibody Subtype/Subclass ELISA Kits

Catalog # 3092, 3093, 3094, 3095, and 3096

For Research Use Only - Not Human or Therapeutic Use

INTRODUCTION

Wheat is the most widely consumed food grain in the world. Wheat proteins are categorized into four fractions based on their solubility in solvents: water (albumins), water containing salt (globulins), alcohol (gliadins), and alkali or acid solution (glutenin) (1). One of these proteins, glutenin is known for its resistance to dissolving in common solvents, as well as for the ability to form aggregates. Glutenin, which can range in MW from 30 to 140 kDa, is divided into two groups: high molecular weight glutenin subunits (HMW: 90–140 kDa), which play a critical role in gluten functionality, and low molecular weight glutenin subunits (LMW: 30–75 kDa), which are partially soluble in 70% ethanol and share similarities with some gliadins (2). Glutenin plays a critical role in activating both the innate and adaptive immune responses, which results in the immune-mediated injury of the intestine such as high intestinal permeability and lamina propria inflammatory cell infiltration. Therefore, glutenin can trigger Celiac Disease (CD) which is an autoimmune disorder in genetically susceptible individuals. Many CD patients even have antibodies which recognize amino acid sequences (QQPGQ, QQPGQQQ, and QQSGQGQ) in HMW-glutenin (3).

Mouse CD models have been widely used to study the pathogenesis of glutenin and the resulting immune responses. The mice who receive glutenin have significantly higher serum glutenin specific IgE and IgG1 antibodies than controls (4, 5). To evaluate the humoral immunity against glutenin in the mouse CD models, Chondrex, Inc. provides ELISA kits for assaying mouse anti-glutenin subtype and subclass antibodies, including IgG, IgG1, IgG2a, IgG2, and IgM antibodies. Chondrex, Inc. also offers ELISA kits for assaying antibody subtypes/subclasses against gliadin, ovalbumin, house dust mite, and peanut allergens, as well as total immunoglobulin subtypes/subclasses levels. For more information, please visit www.chondrex.com or contact support@chondrex.com.

KIT COMPONENTS

	Item	Quantity	Amount	Storage
Standard	IgG (30921) – 100 ng IgG1 (30931) – 100 ng IgG2a (30941) – 100 ng IgG2b (30951) – 200 ng IgM (30961) – 100 ng	1 vial	Lyophilized	-20°C
Secondary Antibody (peroxidase-conjugated polyclonal antibodies)	IgG (30113) IgG1 (30133) IgG2a (30153) IgG2b (203013) IgM (30173)	2 vials	50 µl	-20°C
Solution C - Sample/Standard/Secondary Antibody Dilution Buffer (30106)		2 bottles	50 ml	-20°C
TMB Solution (90023)		2 vials	0.2 ml	-20°C
Chromogen Dilution Buffer (90022)		1 bottle	20 ml	-20°C
Stop Solution - 2N Sulfuric Acid (9016)		1 bottle	10 ml	-20°C
Wash Buffer, 20X (9005)		1 bottle	50 ml	-20°C
Glutenin Coated ELISA Plate (Orange)		1 each	96-well (8-well strips x 12)	-20°C

NOTES BEFORE USING ASSAY

NOTE 1: It is recommended that the standard and samples be run in duplicate.

NOTE 2: Warm up all buffers to room temperature before use.

NOTE 3: Crystals may form in Wash Buffer, 20X when stored at cold temperatures. If crystals have formed, warm the wash buffer by placing the bottle in warm water until crystals are completely dissolved.

NOTE 4: Measure exact volume of buffers using a serological pipet, as extra buffer is provided.

NOTE 5: Cover the plate with plastic wrap or a plate sealer after each step to prevent evaporation from the outside wells of the plate.

NOTE 6: For partial reagent use, please see the assay protocol's corresponding step for the appropriate dilution ratio. For example, if the protocol dilutes 50 μ l of a stock solution in 10 ml of buffer for 12 strips, then for 6 strips, dilute 25 μ l of the stock solution in 5 ml of buffer. Partially used stock reagents may be kept in their original vials and stored at -20°C for use in a future assay.

NOTE 7: This kit contains animal components from non-infectious animals and should be treated as potential biohazards in use and for disposal.

NOTE 8: Depending on the isotypes, subtypes, and targeting epitopes of antibodies, the binding affinity of individual antibodies varies significantly. Therefore, the total IgG antibody concentration calculated as the sum of individual IgG subtypes might not perfectly match the total IgG concentration as determined by the Mouse Anti-Glutenin IgG Antibody ELISA Kit (Cat # 3092).

ASSAY OUTLINE



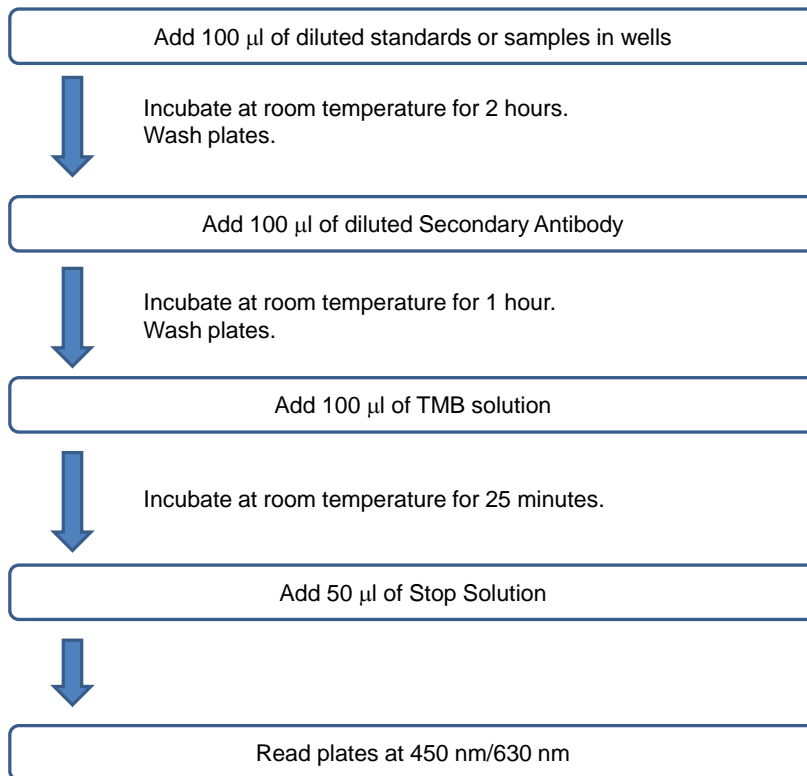


PLATE MAPPING

Example of the Mouse Anti-Glutelin IgG, IgG1, IgG2a, and IgM Antibody ELISA Kits

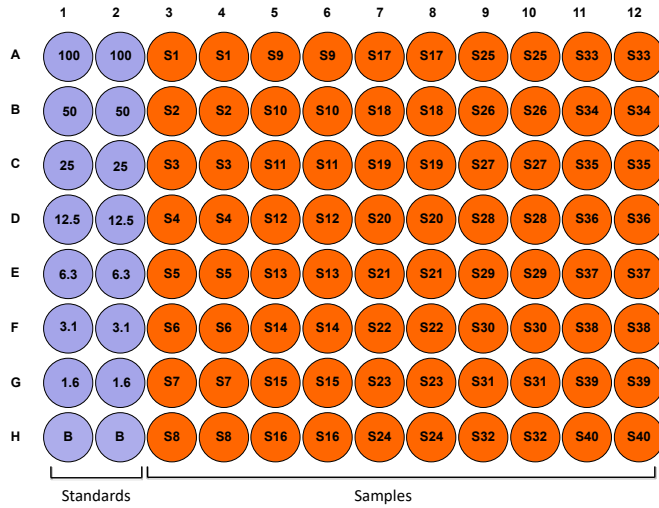


Example of the Mouse Anti-Glutelin IgG2b Antibody ELISA Kits

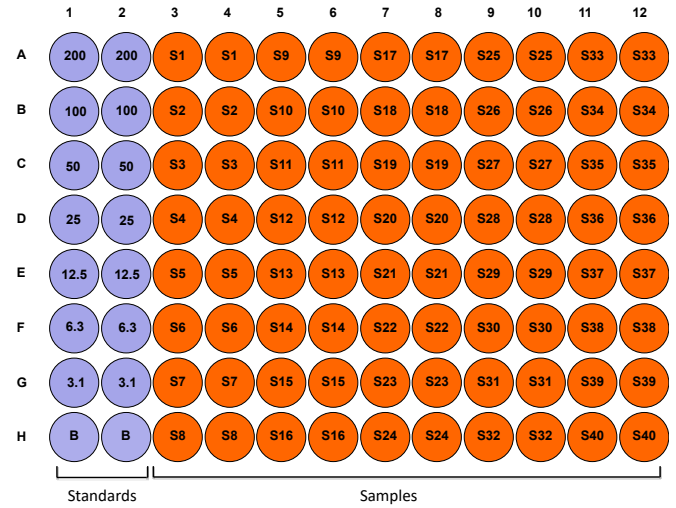

ASSAY PROCEDURE

- Prepare Standard Dilutions:** Please see the figure below for each assay's recommended standard range. Dissolve one vial of Standard in 1 ml of Sample/Standard/Secondary Antibody Dilution Buffer (Solution C) and keep it as a standard stock. Dilute the standard stock accordingly to get the first stock solution. Then serially dilute it with Solution C. For example, mix 250 μ l of the first stock solution with an equal volume of Solution C to make the second stock solution, and then repeat it five more times. The original standard stock solution may be stored at -20°C for use in a future assay. Chondrex, Inc. recommends making fresh serial dilutions for each assay.

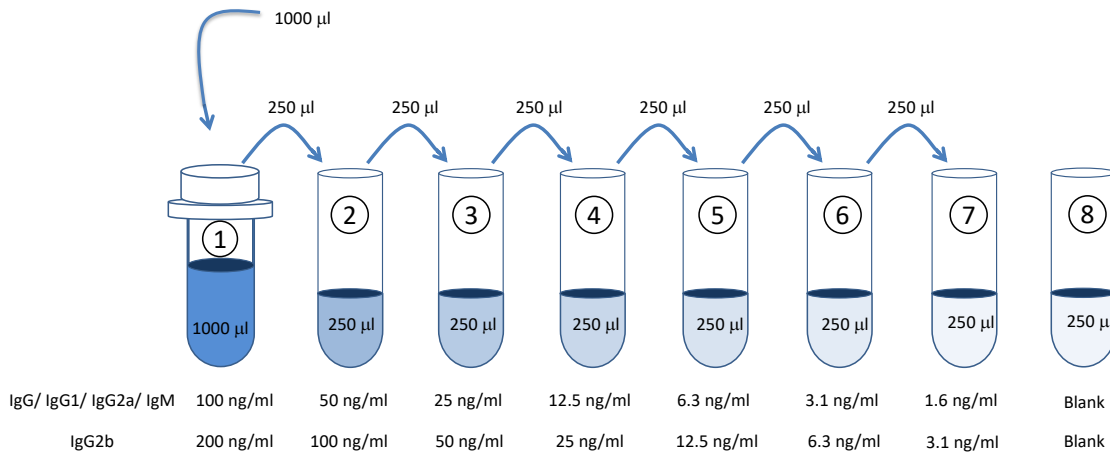

- Prepare Sample Dilutions:** The dilution of serum from mice immunized with glutenin varies (1:1000 or more) depending on the immunization schedule and timing of serum collection. In general, no antibodies against glutenin are observed in normal serum at a 1:1000 dilution.

NOTE: Because regular mouse chows may contain gluten, unimmunized naive mice may have anti-glutenin antibodies in serum.

- Add Standards and Samples:** Add 100 μ l of standards, Solution C (blank), and samples to wells in duplicate. Incubate at room temperature for 2 hours.
- Wash:** Dilute 50 ml of 20X wash buffer in 950 ml of distilled water (1X wash buffer). Wash the plate with 1X wash buffer at least 3 times using a wash bottle with manifold or an automated plate washer. Empty the plate by inverting it and blotting on a paper towel to remove excess liquid. *Do not allow the plate to dry out.*
- Add Secondary Antibody:** Dilute one vial of Secondary Antibody in 10 ml Sample/Standard/Secondary Antibody Dilution Buffer (Solution C). Add 100 μ l of secondary antibody solution to each well and incubate at room temperature for 1 hour.

Strip #	2 nd Antibody (μ l)	Solution C (ml)
2	8	1.7
4	17	3.3
6	25	5.0
8	33	6.6
10	42	8.2
12	50	10.0

- Wash:** Wash the plate with 1X wash buffer at least 3 times using a wash bottle with manifold or an automated plate washer. Empty the plate by inverting it and blotting on a paper towel to remove excess liquid. *Do not allow the plate to dry out.*
- Add TMB Solution:** Use new tubes when preparing TMB solution. Just prior to use, prepare TMB solution with Chromogen Dilution Buffer as shown in the following table. Add 100 μ l of TMB solution to each well immediately after washing the plate and incubate for 25 minutes at room temperature


Strip #	TMB (μ l)	Chromogen Dilution Buffer (ml)
2	34	1.7
4	66	3.3
6	100	5.0
8	132	6.6
10	164	8.2
12	200	10.0

- Stop:** Stop the reaction with 50 μ l of 2N Sulfuric Acid (Stop Solution) to each well.
- Read Plate:** Read the OD values at 450 nm. If the OD values of samples are greater than the OD values of the highest standard, re-assay the samples at a higher dilution. A 630 nm filter can be used as a reference.

CALCULATING RESULTS

- Average the duplicate OD values for the standards, blanks (B), and test samples.
- Subtract the averaged blank OD values from the averaged OD values of the standards and test samples.
- Plot the OD values of the standards against the ng/ml of antibody standard. Using a log/log plot will linearize the data. Figure 1 shows an example of a standard curve for anti-glutenin Ig antibodies.
- The ng/ml of antibody in test samples can be calculated using regression analysis. Multiply it by the sample dilution factor to obtain the antibody concentration (ng/ml) in original test samples.

Figure 1 - Typical Standard Curves for the Anti-Glutenin Antibody ELISA Kits

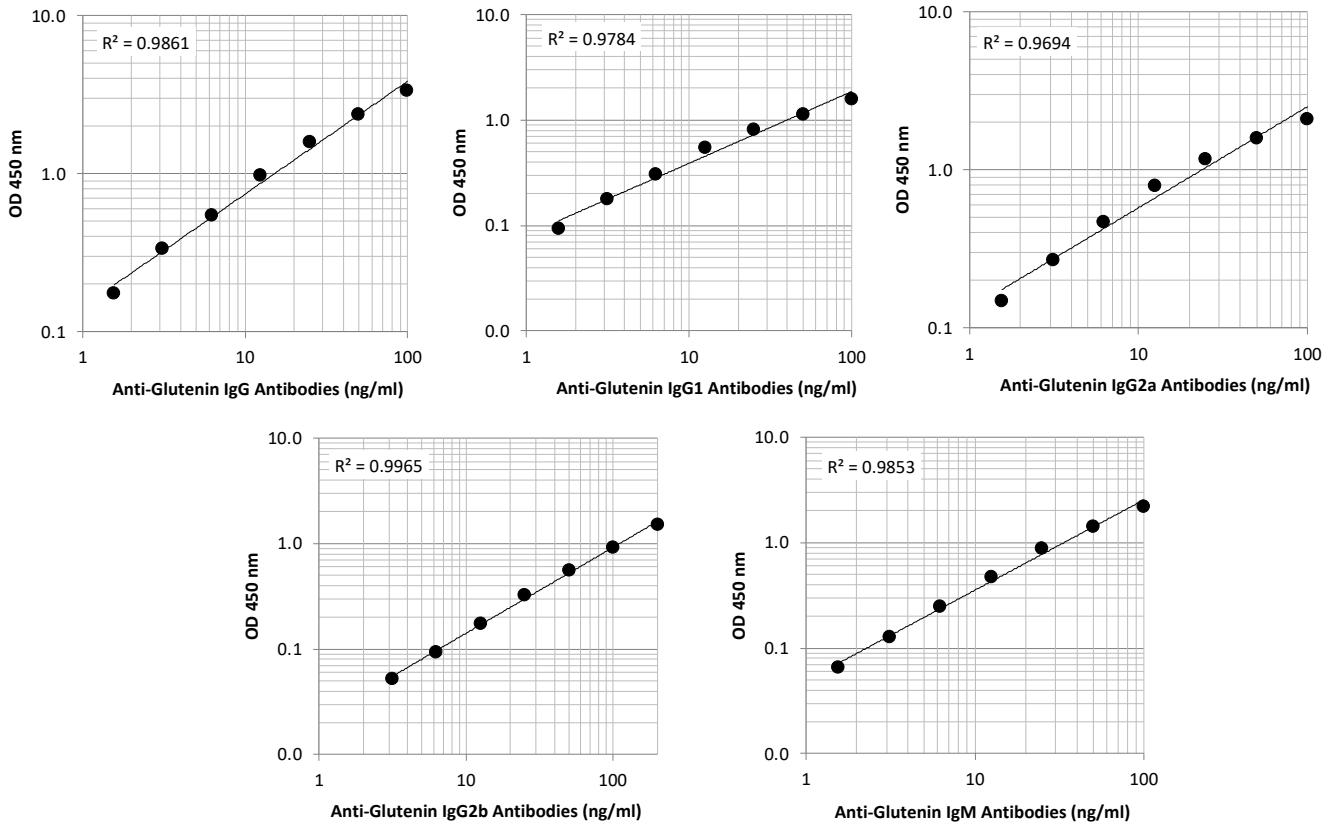

VALIDATION DATA

Table 1 - Reproducibility Data for the Mouse Anti-Glutenin IgG Antibody ELISA Kit

Test	3.1 ng/ml	12.5 ng/ml	50 ng/ml
Intra-Assay CV (%)	3.1	2.6	2.8
Inter-Assay CV (%)	5.3	3.4	4.1
Spike Test* (%)	97%	108%	108%

Table 2 - Reproducibility Data for the Mouse Anti-Glutenin IgG1 Antibody ELISA Kit

Test	3.1 ng/ml	12.5 ng/ml	50 ng/ml
Intra-Assay CV (%)	2.0	3.5	0.9
Inter-Assay CV (%)	10.0	3.7	4.4
Spike Test* (%)	108%	102%	104%

Table 3 - Reproducibility Data for the Mouse Anti-Glutenin IgG2a Antibody ELISA Kit

Test	3.1 ng/ml	12.5 ng/ml	50 ng/ml
Intra-Assay CV (%)	9.1	6.6	6.5
Inter-Assay CV (%)	5.5	1.9	5.9
Spike Test* (%)	103%	102%	98%

Table 4 - Reproducibility Data for the Mouse Anti-Glutenin IgG2b Antibody ELISA Kit

Test	6.3 ng/ml	25 ng/ml	100 ng/ml
Intra-Assay CV (%)	9.4	4.8	8.8
Inter-Assay CV (%)	4.5	5.3	8.4
Spike Test* (%)	103%	105%	97%

Table 5 - Reproducibility Data for the Mouse Anti-Glutenin IgM Antibody ELISA Kit

Test	3.1 ng/ml	12.5 ng/ml	50 ng/ml
Intra-Assay CV (%)	2.3	5.4	7.6
Inter-Assay CV (%)	0.5	5.6	5.3
Spike Test* (%)	99%	98%	109%

*Known amounts of anti-glutenin antibodies were added to samples and then diluted with Sample/Standard/Secondary Antibody Dilution Buffer to assay anti-glutenin antibodies by ELISA.

TROUBLESHOOTING

For frequently asked questions about assays and ELISAs, please see Chondrex, Inc.'s [ELISA FAQ](#) for more information.

REFERENCES

1. N. Inomata, Wheat allergy. *Curr. Opin. Allergy Clin. Immunol.* **9**, 238–243 (2009).
2. H. Wieser, Chemistry of gluten proteins. *Food Microbiol.* **24**, 115–119 (2007).
3. W. Vader, Y. Kooy, V. Van, R. De, D. Harris, W. Benckhuijsen, *et al.*, The gluten response in children with celiac disease is directed toward multiple gliadin and glutenin peptides. *Gastroenterology* **122**(7), 1729-37 (2002).
4. Y. Wang, X. Li, S. Wu, L. Dong, Y. Hu, J. Wang, Y. Zhang, S. Wang, Methylglyoxal decoration of glutenin during heat processing could alleviate the resulting allergic reaction in mice. *Nutrients* **12**, 2844 (2020).
5. H. Kozai, H. Yano, T. Matsuda, Y. Kato, Wheat-dependent exercise-induced anaphylaxis in mice is caused by gliadin and glutenin treatments. *Immunol. Lett.* **102**, 83–90 (2006).